Procedure for Generating a Zone PID Table

Depending on your needs, you can create your own Zone PID table with ten temperature zones.

NOTE: The controller must be in zone mode. This is accomplished by pressing **Control Setup** and then **Next Setting** until you reach "Control Mode". Use ▲ or ▼to select the "zone" option. Press **Save Screen** until you return to the temperature display.

PROCEDURE TO ENTER YOUR OWN TABLE OF PID VALUES:

1) Press **Zone Settings**

You should see the following screen:

Loop: 1 ZONE SETTINGS					
Top 1 100 2 100 3 100 4 100 5 100	P 0:0 0:0 0:0 0:0	I 0.0 0.0 0.0 0.0	000000	Mout % 0.00 0.00 0.00 0.00	Range OFF OFF OFF OFF

- 2) Press **Enter** or **Next Setting** to advance through the entries
- 3) Use the keypad to type in the desired values and push **Enter** to move along
- 4) When you are done, press **Save Screen**

